

Spectroquant®

Prove

The following methods with the corresponding method numbers are programmed into the photometer. The measurements can be made without any further adjustments. In addition several applications (AppNotes) are pre-programmed. The full description of the AppNotes and their procedure you can find on: www.emdmillipore.com/aaf. The applications are grouped into the following segments.

Legend:

- Brewery AppNotes
- Water and Waste Water AppNotes
- ICUMSA and Oil AppNotes
- Further AppNotes



Overview of preprogrammed Methods and AppNotes

| Meth. No. | Determination | Art. No. | Total range | Cell size [mm] | Method |
|--|--|----------|---------------------------------------|----------------|--------------------------------|
| 208 | Acid Capacity Cell Test to pH 4.3 (total alkalinity) | 101758 | 0.40 – 8.00 mmol/l | – | Indicator reaction |
| 2518 | ADMI | AppNote | 2.0 – 100.0 | | Inherent color |
| 2517 | ADMI | AppNote | 10 – 500 | | Inherent color |
| 2612 | α Acids ²⁾ | AppNote | 0 – 80 mg/l | | Inherent color |
| 196 | Aluminium Cell Test ¹⁾ | 100594 | 0.02 – 0.50 mg/l Al | – | Chromazurole S |
| 43 | Aluminium Test ¹⁾ | 114825 | 0.020 – 1.20 mg/l Al | 10, 20, 50 | Chromazurole S |
| Amino nitrogen, free – see Free Amino Nitrogen | | | | | |
| 2520 | Ammonia, free | AppNote | 0.00 – 3.65 mg/l NH ₃ | 10, 20, 50 | as ammonium |
| 104 | Ammonium Cell Test | 114739 | 0.010 – 2.000 mg/l NH ₄ -N | – | Indophenol blue |
| 51 | Ammonium Cell Test | 114558 | 0.20 – 8.00 mg/l NH ₄ -N | – | Indophenol blue |
| 52 | Ammonium Cell Test | 114544 | 0.5 – 16.0 mg/l NH ₄ -N | – | Indophenol blue |
| 53 | Ammonium Cell Test | 114559 | 4.0 – 80.0 mg/l NH ₄ -N | – | Indophenol blue |
| 54 | Ammonium Test | 114752 | 0.010 – 3.00 mg/l NH ₄ -N | 10, 20, 50 | Indophenol blue |
| 155 | Ammonium Test | 100683 | 2.0 – 75.0 mg/l NH ₄ -N | 10 | Indophenol blue |
| 163 | Ammonium Test | 100683 | 5 – 150 mg/l NH ₄ -N | 10 | Indophenol blue |
| 2601 | Anthocyanogenes ²⁾ | AppNote | 0 – 100 mg/l | | Acidic hydrolysis |
| 130 | Antimony in water and wastewater | AppNote | 0.10 – 8.00 mg/l Sb | | Brilliant green |
| 156 | AOX Cell Test ¹⁾ | 100675 | 0.05 – 2.50 mg/l AOX | – | Oxidation to chloride |
| 132 | Arsenic Test ¹⁾ | 101747 | 0.001 – 0.100 mg/l As | 10, 20 | Ag-DDTC |
| 2603 | Bitterness – beer ²⁾ | AppNote | 1 – 80 BU | | UV absorption |
| 2604 | Bitterness – wort ²⁾ | AppNote | 1 – 120 BU | | UV absorption |
| 157 | BOD Cell Test ¹⁾ | 100687 | 0.5 – 3,000 mg/l BOD | – | Modification of Winkler method |
| 164 | Boron Cell Test ¹⁾ | 100826 | 0.05 – 2.00 mg/l B | – | Azomethine H |
| 46 | Boron Test ¹⁾ | 114839 | 0.050 – 0.800 mg/l B | 10 | Rosocyanine |

¹⁾ turbidity correction possible | ²⁾ the analytical procedure for this method is given in the manual of the "Brewery Methods Prove"

³⁾ individual calibration necessary | ⁴⁾ 100 mm cuvette use only for SQ Prove 600

⁵⁾ incapable measurement with SQ Prove 100 | ⁶⁾ difference in test procedure, see package insert



Overview of preprogrammed Methods and AppNotes

| Meth. No. | Determination | Art. No. | Total range | Cell size [mm] | Method |
|--|--|---|--|----------------|--|
| 195 | Bromate in water and drinking water | AppNote | 0.003 – 0.120 mg/l BrO ₃ | – | 3,3'-Dimethylnaphtidine |
| 146 | Bromine Test ¹⁾ | 100605 | 0.020 – 10.00 mg/l Br ₂ | 10 | S-DPD |
| 157 | BSB Cell Test ¹⁾ | 100687 | 0.5 – 3,000 mg/l BSB | – | Modification of Winkler method |
| 67 | Cadmium Cell Test | 114834 | 0.025 – 1.000 mg/l Cd | – | Cadion derivate |
| 183 | Cadmium Test | 101745 | 0.0020 – 0.500 mg/l Cd | 10, 20, 50 | Cadion derivate |
| 165 | Calcium Cell Test ¹⁾ | 100858 | 10 – 250 mg/l Ca | – | Phthalein purple |
| 42 | Calcium Test ^{1), 6)} | 114815 | 5 – 160 mg/l Ca | 10, 20 | Glyoxal-bis-hydroxyanil |
| 125 | Calcium Test sensitive ^{1), 6)} | 114815 | 1.0 – 15.0 mg/l Ca | 10 | Glyoxal-bis-hydroxyanil |
| 304 | Calcium Test ³⁾ | 100049 | 0.20 – 4.00 mg/l Ca | 10 | Phthalein derivate |
| Carbohydrates, total – see Total Carbohydrates | | | | | |
| 2523 | Carotene (palm oil) | AppNote | 10 – 7,500 mg/kg | – | Inherent color |
| 95 | Chloride Cell Test ¹⁾ | 114730 | 5 – 125 mg/l Cl | – | Iron(III)-thiocyanat |
| 110 | Chloride Test ^{1), 6)} | 114897 | 2.5 – 25.0 mg/l Cl | 10 | Iron(III)-thiocyanat |
| 63 | Chloride Test ^{1), 6)} | 114897 | 10 – 250 mg/l Cl | 10 | Iron(III)-thiocyanat |
| 218 | Chloride Cell Test ¹⁾ | 101804 | 0.5 – 15.0 mg/l Cl | – | Iron(III)-thiocyanat |
| 219 | Chloride Test ¹⁾ | 101807 | 0.10 – 5.00 mg/l Cl | 50 | Iron(III)-thiocyanat |
| 141 | Chlorine Cell Test ¹⁾ (free chlorine) | 100595 | 0.03 – 6.00 mg/l Cl ₂ | – | S-DPD |
| 142 | Chlorine Cell Test ¹⁾ (free chlorine + total chlorine) | 100597 | 0.03 – 6.00 mg/l Cl ₂ | – | S-DPD |
| 143 | Chlorine Test ¹⁾ (free chlorine) | 100598 | 0.010 – 6.00 mg/l Cl ₂ | 10, 20, 50 | S-DPD |
| 145 | Chlorine Test ¹⁾ (total chlorine) | 100602 | 0.010 – 6.00 mg/l Cl ₂ | 10, 20, 50 | S-DPD |
| 144 | Chlorine Test ¹⁾ (free chlorine + total chlorine) | 100599 | 0.010 – 6.00 mg/l Cl ₂ | 10, 20, 50 | S-DPD |
| 194 | Chlorine Cell Test ¹⁾ (free chlorine + total chlorine) | 100086/ 100087/ 100088/ 100089 | 0.03 – 6.00 mg/l Cl ₂ | 10 | DPD |
| 306 | Chlorine Test ¹⁾ (free chlorine + total chlorine) | 100086/ 100087/ 100088 | 0.010 – 1.000 mg/l Cl ₂ | 10 | DPD |
| 149 | Chlorine Dioxide Test ¹⁾ | 100608 | 0.020 – 10.00 mg/l ClO ₂ | 10 | S-DPD |
| 2509 | Chlorophyll-a (DIN/ISO) | AppNote | result in µg/l Chl-a or Phaeo | – | Inherent color |
| 2504 | Chlorophyll-a (APHA/ASTM) | AppNote | result in mg/m ³ Chl-a or Phaeo | – | Inherent color |
| 2507 | Chlorophyll-a, -b, -c (APHA/ASTM) | AppNote | result in mg/m ³ Chl-a, -b, -c | – | Inherent color |
| 39 | Chromate Cell Test ¹⁾ | 114552 | 0.05 – 2.00 mg/l Cr | – | Diphenylcarbazide |
| 39 | Chromate Cell Test ¹⁾ (total chromium) | 114552 | 0.05 – 2.00 mg/l Cr | – | Peroxodisulfate oxidation / Diphenylcarbazide |
| 40 | Chromate Test ¹⁾ | 114758 | 0.010 – 3.00 mg/l Cr | 10, 20, 50 | Diphenylcarbazide |
| 20 | Chromium Baths | AppNote | 4.0 – 400 g/l CrO ₃ | – | Inherent color |
| 305 | Cobalt in water | AppNote | 0.5 – 10.0 mg/l Co | – | Nitroso-R salt |
| 31 | COD Cell Test ¹⁾ | 114560 | 4.0 – 40.0 mg/l COD | – | Chromosulfuric acid oxidation / chromate determination |
| 211 | COD Cell Test ¹⁾ | 101796 | 5.0 – 80.0 mg/l COD | – | Chromosulfuric acid oxidation / chromate determination |

¹⁾ turbidity correction possible | ²⁾ the analytical procedure for this method is given in the manual of the "Brewery Methods Prove"

³⁾ individual calibration necessary | ⁴⁾ 100 mm cuvette use only for SQ Prove 600

⁵⁾ incapable measurement with SQ Prove 100 | ⁶⁾ difference in test procedure, see pagacke insert



Overview of preprogrammed Methods and AppNotes

| Meth. No. | Determination | Art. No. | Total range | Cell size [mm] | Method |
|-----------|---|----------|------------------------------|----------------|--|
| 14 | COD Cell Test ¹⁾ | 114540 | 10 – 150 mg/l COD | – | Chromosulfuric acid oxidation / chromate determination |
| 105 | COD Cell Test ¹⁾ | 114895 | 15 – 300 mg/l COD | – | Chromosulfuric acid oxidation / chromate determination |
| 93 | COD Cell Test ¹⁾ | 114690 | 50 – 500 mg/l COD | – | Chromosulfuric acid oxidation / chromate determination |
| 23 | COD Cell Test ¹⁾ | 114541 | 25 – 1,500 mg/l COD | – | Chromosulfuric acid oxidation / chromium(III) determination |
| 94 | COD Cell Test ¹⁾ | 114691 | 300 – 3,500 mg/l COD | – | Chromosulfuric acid oxidation / chromium(III) determination |
| 24 | COD Cell Test ¹⁾ | 114555 | 500 – 10,000 mg/l COD | – | Chromosulfuric acid oxidation / chromium(III) determination |
| 209 | COD Cell Test ¹⁾ | 101797 | 5,000 – 90,000 mg/l COD | – | Chromosulfuric acid oxidation / chromium(III) determination |
| 137 | COD Cell Test (Hg free) ¹⁾ | 109772 | 10 – 150 mg/l COD | – | Chromosulfuric acid oxidation / chromate determination |
| 138 | COD Cell Test (Hg free) ¹⁾ | 109773 | 100 – 1,500 mg/l COD | – | Chromosulfuric acid oxidation / chromium(III) determination |
| 220 | COD Cell Test for seawater ¹⁾ | 117058 | 5.0 – 60.0 mg/l COD | – | Chloride depletion / chromosulfuric acid oxidation / chromate determination |
| 221 | COD Cell Test for seawater ¹⁾ | 117059 | 50 – 3,000 mg/l COD | – | Chloride depletion / chromosulfuric acid oxidation / chromium(III) determination |
| 15 | Color α (436) (spectral absorptions coefficient) | AppNote | 0.1 – 250 m ⁻¹ | | Measurement at 436 nm |
| 61 | Color α (525) (spectral absorptions coefficient) | AppNote | 0.1 – 250 m ⁻¹ | | Measurement at 525 nm |
| 78 | Color α (620) (spectral absorptions coefficient) | AppNote | 0.1 – 250 m ⁻¹ | | Measurement at 620 nm |
| 303 | Color (410) (EN 7887) | AppNote | 2 – 2,500 mg/l Pt | | Measurement at 410 nm |
| 2602 | Color – EBC ²⁾ | AppNote | 0.0 – 60.0 EBC Units | | Inherent color |
| 32 | Color Hazen ¹⁾ | AppNote | 0.2 – 500 mg/l Pt/Co (Hazen) | | Platinum-cobalt-Standard Method, measurement at 340 nm |
| 179 | Color Hazen ¹⁾ | AppNote | 0 – 1,000 mg/l Pt/Co (Hazen) | | Platinum-cobalt-Standard Method, measurement at 445 nm |
| 180 | Color Hazen ¹⁾ | AppNote | 0 – 1,000 mg/l Pt/Co (Hazen) | | Platinum-cobalt-Standard Method, measurement at 455 nm |
| 181 | Color Hazen ¹⁾ | AppNote | 0 – 1,000 mg/l Pt/Co (Hazen) | | Platinum-cobalt-Standard Method, measurement at 465 nm |

¹⁾ turbidity correction possible | ²⁾ the analytical procedure for this method is given in the manual of the "Brewery Methods Prove"

³⁾ individual calibration necessary | ⁴⁾ 100 mm cuvette use only for SQ Prove 600

⁵⁾ incapable measurement with SQ Prove 100 | ⁶⁾ difference in test procedure, see package insert



Overview of preprogrammed Methods and AppNotes

| Meth. No. | Determination | Art. No. | Total range | Cell size [mm] | Method |
|---|--|----------|---|----------------|---|
| Color of sugar solutions – see ICUMSA Color | | | | | |
| 2613 | Copper – EBC ²⁾ | AppNote | 0.10 – 5.00 mg/l Cu | | Cuprethol |
| 26 | Copper Cell Test ¹⁾ | 114553 | 0.05 – 8.00 mg/l Cu | – | Cuprizone |
| 27 | Copper Test ¹⁾ | 114767 | 0.02 – 6.00 mg/l Cu | 10, 20, 50 | Cuprizone |
| 83 | Copper Baths | AppNote | 2.0 – 80.0 g/l Cu | | Inherent color |
| 228 | Cyanide Cell Test ¹⁾ (free cyanide) | 102531 | 0.010 – 0.500 mg/l CN | – | Barbituric acid + pyridinecarboxylic acid |
| 75 | Cyanide Cell Test ¹⁾ (free cyanide) | 114561 | 0.010 – 0.500 mg/l CN | – | Barbituric acid + pyridinecarboxylic acid |
| 75 | Cyanide Cell Test ¹⁾ (readily liberated cyanide) | 114561 | 0.010 – 0.500 mg/l CN | – | Citric acid / barbituric acid + pyridinecarboxylic acid |
| 109 | Cyanide Test ¹⁾ (free cyanide) | 109701 | 0.0020 – 0.500 mg/l CN | 10, 20, 50 | Barbituric acid + pyridinecarboxylic acid |
| 109 | Cyanide Test ¹⁾ (readily liberated cyanide) | 109701 | 0.0020 – 0.500 mg/l CN | 10, 20, 50 | Citric acid / barbituric acid + pyridinecarboxylic acid |
| 210 | Cyanuric Acid Test | 119253 | 2 – 160 mg/l Cyan Acid | 20 | Triazine derivative |
| 2528 | delta K268 (olive oil) | AppNote | -0.10 – 1.00 | | UV absorption |
| 2529 | delta K270 (olive oil) | AppNote | -0.10 – 1.00 | | UV absorption |
| Diacetyl – see Vicinal Diketones | | | | | |
| 2524 | DOBI (palm oil) | AppNote | 0.00 – 4.00 | | UV absorption |
| 2626 | Flavonoids ²⁾ | AppNote | 3.0 – 200.0 mg/l | | 4-Dimethylaminocinnamaldehyde |
| 215 | Fluoride Cell Test ¹⁾ | 100809 | 0.10 – 1.80 mg/l F | – | Alizarin complexone |
| 216 | Fluoride Cell Test sensitive | 100809 | 0.025 – 0.500 mg/l F | 50 | Alizarin complexone |
| 166 | Fluoride Test ^{1), 6)} | 114598 | 0.10 – 2.00 mg/l F | 10 | Alizarin complexone |
| 167 | Fluoride Test ^{1), 6)} | 114598 | 1.0 – 20.0 mg/l F | 10 | Alizarin complexone |
| 217 | Fluoride Test | 100822 | 0.02 – 2.00 mg/l F | 50 | SPADNS |
| 28 | Formaldehyde Cell Test ¹⁾ | 114500 | 0.10 – 8.00 mg/l HCHO | – | Chromotropic acid |
| 91 | Formaldehyde Test ¹⁾ | 114678 | 0.02 – 8.00 mg/l HCHO | 10, 20, 50 | Chromotropic acid |
| 2606 | Free Amino Nitrogen beer / wort ²⁾ | AppNote | 0 – 400 mg/l | | Ninhydrin |
| 45 | Gold Test | 114821 | 0.5 – 12.0 mg/l Au | 10 | Rhodamine B |
| Hardness – see Total Hardness or Residual Hardness | | | | | |
| Hazen – see Color Hazen | | | | | |
| 44 | Hydrazine Test ¹⁾ | 109711 | 0.005 – 2.00 mg/l N ₂ H ₄ | 10, 20, 50 | 4-Dimethylaminobenzaldehyde |
| 99 | Hydrogen Peroxide Cell Test ¹⁾ | 114731 | 2.0 – 20.0 mg/l H ₂ O ₂ | – | Titanyl sulfate |
| 128 | Hydrogen Peroxide Cell Test sensitive ¹⁾ | 114731 | 0.25 – 5.00 mg/l H ₂ O ₂ | 50 | Titanyl sulfate |
| 198 | Hydrogen Peroxide Test | 118789 | 0.015 – 6.00 mg/l H ₂ O ₂ | 10, 20 | Phenanthroline derivative |
| 2548 | ICUMSA Color GS1/3-7 | AppNote | 0 – 50,000 IU _{7,0} | | Inherent color |
| 2549 | ICUMSA Color GS2/3-9 | AppNote | 0 – 600 IU _{7,0} | | Inherent color |
| 2550 | ICUMSA Color GS2/3-10 | AppNote | 0 – 50 IU _{7,0} | | Inherent color |
| 2551 | ICUMSA Color GS9/1/2/3-8 | AppNote | 0 – 20,000 IU _{7,0} | | Inherent color |
| 147 | Iodine Test ¹⁾ | 100606 | 0.050 – 10.00 mg/l I ₂ | 10, 20, 50 | S-DPD |
| 2615 | Iodine Test, photometric ²⁾ | AppNote | 0.00 – 0.80 | | Iodine |
| 2616 | Iodine Test, photometric ²⁾ | AppNote | 0.00 – 0.80 | | Iodine |
| 33 | Iodine Color Number | AppNote | 0.010 – 3.00 | | Measurement at 340 nm |
| 21 | Iodine Color Number | AppNote | 0.2 – 50.0 | | Measurement at 445 nm |
| 2623 | Iron – EBC ²⁾ | AppNote | 0.000 – 1.000 mg/l Fe | | Triazine |
| 2624 | Iron – EBC ²⁾ | AppNote | 0.000 – 0.800 mg/l Fe | | Triazine |
| 37 | Iron Cell Test | 114549 | 0.05 – 4.00 mg/l Fe | – | Triazine |

¹⁾ turbidity correction possible | ²⁾ the analytical procedure for this method is given in the manual of the "Brewery Methods Prove"

³⁾ individual calibration necessary | ⁴⁾ 100 mm cuvette use only for SQ Prove 600

⁵⁾ incapable measurement with SQ Prove 100 | ⁶⁾ difference in test procedure, see package insert



Overview of preprogrammed Methods and AppNotes

| Meth. No. | Determination | Art. No. | Total range | Cell size [mm] | Method |
|--|---|----------|---|-------------------------------|--|
| 106 | Iron Cell Test ¹⁾ | 114896 | 1.0 – 50.0 mg/l Fe (Fe(II) and Fe(III)) | | 2,2'-Dipyridyl |
| 38 | Iron Test | 114761 | 0.005 – 5.00 mg/l Fe | 10, 20, 50, 100 ⁴⁾ | Triazine |
| 161 | Iron Test ¹⁾ | 100796 | 0.010 – 5.00 mg/l Fe (Fe(II) and Fe(III)) | 10, 20, 50 | 1,10-Phenanthroline |
| 2611 | Iso- α Acids ²⁾ | AppNote | 0 – 60 | | UV absorption |
| 2525 | K232 (olive oil) | AppNote | 0.00 – 4.00 | | UV absorption |
| 2526 | K268 (olive oil) | AppNote | 0.00 – 4.00 | | UV absorption |
| 2527 | K270 (olive oil) | AppNote | 0.00 – 4.00 | | UV absorption |
| 66 | Lead Cell Test ¹⁾ | 114833 | 0.10 – 5.00 mg/l Pb | – | PAR |
| 160 | Lead Test ¹⁾ | 109717 | 0.010 – 5.00 mg/l Pb | 10, 20, 50 | PAR |
| 158 | Magnesium Cell Test ¹⁾ | 100815 | 5.0 – 75.0 mg/l Mg | – | Phthalein purple |
| 159 | Manganese Cell Test ¹⁾ | 100816 | 0.10 – 5.00 mg/l Mn | – | Formaldehyde |
| 19 | Manganese Test ¹⁾ | 114770 | 0.010 – 10.00 mg/l Mn | 10, 20, 50 | Formaldehyde |
| 226 | Manganese Test ¹⁾ | 101846 | 0.005 – 2.00 mg/l Mn | 10, 20, 50 | PAN |
| 135 | Mercury in water and wastewater | AppNote | 0.025 – 1.000 mg/l Hg | | Michler's ketone |
| 175 | Molybdenum Cell Test | 100860 | 0.02 – 1.00 mg/l Mo | – | Brompyrogallol red |
| 206 | Molybdenum Test | 119252 | 0.5 – 45.00 mg/l Mo | 20 | Mercaptoacetic acid |
| 185 | Monochloramine Test | 101632 | 0.050 – 10.00 mg/l Cl ₂ | 10, 20, 50 | Indophenol blue |
| 2614 | Nickel – EBC ²⁾ | AppNote | 0.00 – 5.00 mg/l Ni | | Dimethylglyoxime |
| 17 | Nickel Cell Test ¹⁾ | 114554 | 0.10 – 6.00 mg/l Ni | – | Dimethylglyoxime |
| 18 | Nickel Test ¹⁾ | 114785 | 0.02 – 5.00 mg/l Ni | 10, 20, 50 | Dimethylglyoxime |
| 57 | Nickel Bath | AppNote | 2.0 – 120 g/l Ni | | Inherent color |
| 59 | Nitrate Cell Test ¹⁾ | 114542 | 0.5 – 18.0 mg/l NO ₃ -N | – | Nitrospectral |
| 30 | Nitrate Cell Test ¹⁾ | 114563 | 0.5 – 25.0 mg/l NO ₃ -N | – | 2,6-Dimethylphenol |
| 107 | Nitrate Cell Test ¹⁾ | 114764 | 1.0 – 50.0 mg/l NO ₃ -N | – | 2,6-Dimethylphenol |
| 151 | Nitrate Cell Test ¹⁾ | 100614 | 23 – 225 mg/l NO ₃ -N | – | 2,6-Dimethylphenol |
| 60 | Nitrate Test ¹⁾ | 114773 | 0.2 – 20.0 mg/l NO ₃ -N | 10, 20 | Nitrospectral |
| 139 | Nitrate Cell Test ¹⁾ | 109713 | 0.10 – 25.0 mg/l NO ₃ -N | 10, 20, 50 | 2,6-Dimethylphenol |
| 72 | Nitrate Cell Test in seawater ¹⁾ | 114556 | 0.10 – 3.00 mg/l NO ₃ -N | – | Resorcine |
| 140 | Nitrate Test in seawater ¹⁾ | 114942 | 0.2 – 17.0 mg/l NO ₃ -N | 10 | Resorcine |
| 227 | Nitrate Test | 101842 | 0.3 – 30.0 mg/l NO ₃ -N | 50 | Reduction / Benzoic acid derivative |
| 2503 | Nitrate (UV) | AppNote | 0.0 – 7.0 mg/l NO ₃ -N | | Direct measurement in the UV range |
| 35 | Nitrite Cell Test ¹⁾ | 114547 | 0.010 – 0.700 mg/l NO ₂ -N | – | Griess reaction |
| 197 | Nitrite Cell Test ¹⁾ | 100609 | 1.0 – 90.0 mg/l NO ₂ -N | – | Iron(II)-ethylenediammonium sulfate |
| 36 | Nitrite Test ¹⁾ | 114776 | 0.002 – 1.00 mg/l NO ₂ -N | 10, 20, 50 | Griess reaction |
| 68 | Nitrogen (total) Cell Test | 114537 | 0.5 – 15.0 mg/l N | – | Peroxodisulfate oxidation / Nitrospectral |
| 153 | Nitrogen (total) Cell Test | 100613 | 0.5 – 15.0 mg/l N | – | Peroxodisulfate oxidation / 2,6-Dimethylphenol |
| 108 | Nitrogen (total) Cell Test | 114763 | 10 – 150 mg/l N | – | Peroxodisulfate oxidation / 2,6-Dimethylphenol |
| Oils – see K (olive oil), delta K (olive oil), Carotene (palm oil) or DOBI (palm oil) | | | | | |
| 92 | Oxygen Cell Test ¹⁾ | 114694 | 0.5 – 12.0 mg/l O ₂ | – | Modification of Winkler method |
| 207 | Oxygen Scavengers Test | 119251 | 0.020 – 0.500 mg/l DEHA | 20 | FerroZine® |
| 148 | Ozone Test ¹⁾ | 100607 | 0.010 – 4.00 mg/l O ₃ | 10, 20, 50 | S-DPD |
| 133 | Palladium in water and wastewater | AppNote | 0.05 – 1.25 mg/l Pd | | Thio-Michler's ketone |

¹⁾ turbidity correction possible | ²⁾ the analytical procedure for this method is given in the manual of the "Brewery Methods Prove"

³⁾ individual calibration necessary | ⁴⁾ 100 mm cuvette use only for SQ Prove 600

⁵⁾ incapable measurement with SQ Prove 100 | ⁶⁾ difference in test procedure, see package insert



Overview of preprogrammed Methods and AppNotes

| Meth. No. | Determination | Art. No. | Total range | Cell size [mm] | Method |
|---|---|----------|---|-------------------------------|---|
| 2,3-Pentandion – see Vicinal Diketones | | | | | |
| 186 | pH Cell Test | 101744 | 6.4 – 8.8 | – | Phenol red |
| Phaeophytin (DIN/ISO) / (APHA/ASTM) – see Chlorophyll-a (DIN/ISO) or (APHA/ASTM) | | | | | |
| 73 | Phenol Cell Test ¹⁾ | 114551 | 0.10 – 2.50 mg/l C ₆ H ₅ OH | – | MBTH |
| 176 | Phenol Test ^{1), 6)} | 100856 | 0.025 – 5.00 mg/l C ₆ H ₅ OH | 10, 20, 50 | Aminoantipyrine |
| 177 | Phenol Test ^{1), 6)} | 100856 | 0.002 – 0.100 mg/l C ₆ H ₅ OH | 20 | Aminoantipyrine by extraction |
| Phenols, steam-volatile – see steam-volatiles Phenols | | | | | |
| 212 | Phosphate Cell Test | 100474 | 0.05 – 5.00 mg/l PO ₄ -P | – | Phosphormolybdenum blue |
| 55 | Phosphate Cell Test | 114543 | 0.05 – 5.00 mg/l PO ₄ -P | – | Phosphormolybdenum blue |
| 55 | Phosphate Cell Test (total phosphorus) | 114543 | 0.05 – 5.00 mg/l P | – | Peroxodisulfate oxidation / phosphormolybdenum blue |
| 213 | Phosphate Cell Test | 100475 | 0.5 – 25.0 mg/l PO ₄ -P | – | Phosphormolybdenum blue |
| 86 | Phosphate Cell Test | 114729 | 0.5 – 25.0 mg/l PO ₄ -P | – | Phosphormolybdenum blue |
| 86 | Phosphate Cell Test (total phosphorus) | 114729 | 0.5 – 25.0 mg/l P | – | Peroxodisulfate oxidation / phosphormolybdenum blue |
| 152 | Phosphate Cell Test | 100616 | 3.0 – 100.0 mg/l PO ₄ -P | – | Phosphormolybdenum blue |
| 214 | Phosphate Cell Test | 100673 | 3.0 – 100.0 mg/l PO ₄ -P | – | Phosphormolybdenum blue |
| 214 | Phosphate Cell Test (total phosphorus) | 100673 | 3.0 – 100.0 mg/l P | – | Peroxodisulfate oxidation / phosphormolybdenum blue |
| 56 | Phosphate Test | 114848 | 0.010 – 5.00 mg/l PO ₄ -P | 10, 20, 50, 100 ⁴⁾ | Phosphormolybdenum blue |
| 162 | Phosphate Test | 100798 | 1.0 – 100.0 mg/l PO ₄ -P | 10 | Phosphormolybdenum blue |
| 69 | Phosphate Cell Test ¹⁾ | 114546 | 0.5 – 25.0 mg/l PO ₄ -P | – | Vanadatомolybdate |
| 70 | Phosphate Test ¹⁾ | 114842 | 0.5 – 30.0 mg/l PO ₄ -P | 10, 20 | Vanadatомolybdate |
| Photometric iodine test – see Iodine Test, photometric | | | | | |
| 134 | Platinum in water and wastewater | AppNote | 0.10 – 1.25 mg/l Pt | – | o-Phenylendiamine |
| 103 | Potassium Cell Test | 114562 | 5.0 – 50.0 mg/l K | – | Kalignost®, turbidimetric |
| 150 | Potassium Cell Test | 100615 | 30 – 300 mg/l K | – | Kalignost®, turbidimetric |
| 2617 | Reducing Power ²⁾ | AppNote | 0 – 100 % | – | DPI |
| 98 | Residual Hardness Cell Test ¹⁾ | 114683 | 0.50 – 5.00 mg/l Ca | – | Phthalein |
| 79 | Silicate (Silicic acid) Test | 114794 | 0.11 – 10.70 mg/l SiO ₂ | 10, 20 | Silicomolybdenum blue |
| 81 | Silicate (Silicic acid) Test | 114794 | 0.011 – 1.600 mg/l SiO ₂ | 50 | Silicomolybdenum blue |
| 169 | Silicate (Silicic acid) Test ^{1), 6)} | 100857 | 1.1 – 107.0 mg/l SiO ₂ | 10 | Molybdatosilicate |
| 171 | Silicate (Silicic acid) Test ^{1), 6)} | 100857 | 11 – 1070 mg/l SiO ₂ | 10 | Molybdatosilicate |
| 225 | Silicate (Silicic acid) Test | 101813 | 0.0005 – 0.5000 mg/l SiO ₂ | 50, 100 ⁴⁾ | Silicomolybdenum blue |
| 47 | Silver Test ¹⁾ | 114831 | 0.25 – 3.00 mg/l Ag | 10, 20 | Eosine / 1,10-Phenanthroline |
| 168 | Sodium Cell Test in nutrient solutions ¹⁾ | 100885 | 10 – 300 mg/l Na | – | indirectly as chloride |
| 300 | Spectral Absorption ⁵⁾ Coefficient α(254) | AppNote | 0.5 – 250 m ⁻¹ | – | Measurement at 254 nm |
| 302 | Spectral Absorption Coefficient α(436) | AppNote | 0.5 – 250 m ⁻¹ | – | Measurement at 436 nm |
| 301 | Spectral Attenuation Coefficient μ(254) ^{1), 5)} | AppNote | 0.5 – 250 m ⁻¹ | – | Measurement at 254 nm |
| 2621 | Steam-volatiles Phenols – malt ²⁾ | AppNote | 0.00 – 3.00 mg/kg | – | Aminoantipyrine by extraction |
| 2621 | Steam-volatiles Phenols – beer ²⁾ | AppNote | 0.00 – 0.30 mg/kg | – | Aminoantipyrine by extraction |
| 2622 | Steam-volatiles Phenols – malt ²⁾ | AppNote | 0.00 – 3.00 mg/kg | – | Aminoantipyrine by extraction |
| 2622 | Steam-volatiles Phenols – beer ²⁾ | AppNote | 0.00 – 0.30 mg/kg | – | Aminoantipyrine by extraction |

¹⁾ turbidity correction possible | ²⁾ the analytical procedure for this method is given in the manual of the "Brewery Methods Prove"

³⁾ individual calibration necessary | ⁴⁾ 100 mm cuvette use only for SQ Prove 600

⁵⁾ incapable measurement with SQ Prove 100 | ⁶⁾ difference in test procedure, see pagacke insert



Overview of preprogrammed Methods and AppNotes

| Meth. No. | Determination | Art. No. | Total range | Cell size [mm] | Method |
|---|--|----------|---|----------------|---------------------------------------|
| Sugar solutions, Color of – see ICUMSA Color | | | | | |
| 229 | Sulfate Cell Test | 102532 | 1.0 – 50.0 mg/l SO ₄ | – | Bariumsulfate, turbidimetric |
| 64 | Sulfate Cell Test | 114548 | 5 – 250 mg/l SO ₄ | – | Bariumsulfate, turbidimetric |
| 154 | Sulfate Cell Test | 100617 | 50 – 500 mg/l SO ₄ | – | Bariumsulfate, turbidimetric |
| 82 | Sulfate Cell Test | 114564 | 100 – 1,000 mg/l SO ₄ | – | Bariumsulfate, turbidimetric |
| 65 | Sulfate Test ¹⁾ | 114791 | 25 – 300 mg/l SO ₄ | 10 | Tannin |
| 224 | Sulfate Test | 101812 | 0.50 – 50.0 mg/l SO ₄ | 10, 20, 50 | Bariumsulfate, turbidimetric |
| 230 | Sulfate Test | 102537 | 5 – 300 mg/l SO ₄ | – | Bariumsulfate, turbidimetric |
| 80 | Sulfide Test ¹⁾ | 114779 | 0.020 – 1.50 mg/l S | 10, 20, 50 | Dimethyl-p- phenylendiamin |
| 71 | Sulfite Cell Test ¹⁾ | 114394 | 1.0 – 20.0 mg/l SO ₃ | – | Ellman's reagent |
| 127 | Sulfite Cell Test sensitive ¹⁾ | 114394 | 0.05 – 3.00 mg/l SO ₃ | 50 | Ellman's reagent |
| 187 | Sulfite Test ¹⁾ | 101746 | 1.0 – 60.0 mg/l SO ₃ | 10 | Ellman's reagent |
| 231 | Surfactants (anionic) Cell Test | 102552 | 0.05 – 2.00 mg/l MBAS (methylene blue active substances) | – | Methylene blue |
| 192 | Surfactants (cationic) Cell Test ¹⁾ | 101764 | 0.05 – 1.50 mg/l k-Ten | – | Disulfine blue |
| 193 | Surfactants (nonionic) Cell Test ¹⁾ | 101787 | 0.10 – 7.50 mg/l n-Ten | – | TBPE |
| 182 | Suspended Solids | AppNote | 25 – 750 mg/l SusS | – | |
| 2619 | Thiobarbituric Acid Number ²⁾ | AppNote | 0 – 250 | – | Thiobarbituric acid |
| 100 | Tin Cell Test ¹⁾ | 114622 | 0.10 – 2.50 mg/l Sn | – | Pyrocatechol violet |
| 172 | TOC Cell Test | 114878 | 5.0 – 80.0 mg/l TOC | – | Peroxodisulfate oxidation / Indicator |
| 173 | TOC Cell Test | 114879 | 50 – 800 mg/l TOC | – | Peroxodisulfate oxidation / Indicator |
| 2625 | Total Carbohydrates ²⁾ | AppNote | 0.000 – 6.000 g/100 ml | – | Anthrone |
| 178 | Total Hardness Cell Test ¹⁾ | 100961 | 5 – 215 mg/l Ca | – | Phthalein purple |
| 2610 | Total Polyphenols ²⁾ | AppNote | 1 – 800 mg/l | – | Iron(III) |
| 77 | Turbidity | AppNote | 1 – 100 FAU | – | Measurement at 550 nm |
| 2620 | Vicinal Diketones ²⁾ | AppNote | 0.00 – 1.00 mg/kg | – | Phenylendiamin |
| 222 | Volatile Organic Acids Cell Test ¹⁾ | 101749 | 50 – 3,000 mg/l CH ₃ COOH | – | Esterification |
| 223 | Volatile Organic Acids Test ¹⁾ | 101809 | 50 – 3,000 mg/l CH ₃ COOH | – | Esterification |
| Water hardness – see Total Hardness or Residual Hardness | | | | | |
| 174 | Zinc Cell Test | 100861 | 0.025 – 1.000 mg/l Zn | – | PAR |
| 74 | Zinc Cell Test | 114566 | 0.20 – 5.00 mg/l Zn | – | PAR |
| 41 | Zinc Test ¹⁾ | 114832 | 0.05 – 2.50 mg/l Zn | 10 | Cl-PAN |

¹⁾ turbidity correction possible | ²⁾ the analytical procedure for this method is given in the manual of the "Brewery Methods Prove"

³⁾ individual calibration necessary | ⁴⁾ 100 mm cuvette use only for SQ Prove 600

⁵⁾ incapable measurement with SQ Prove 100 | ⁶⁾ difference in test procedure, see pagacke insert



We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose. Spectroquant® is a trademark of Merck KGaA, Darmstadt, Germany.



For further information on EMD Millipore
and our products contact:

EMD Millipore Corporation
290 Concord Road
Billerica, MA 01821, U.S.A.
www.emdmillipore.com/test-kits

© 2015 Merck KGaA, Darmstadt, Germany. All rights reserved.



www.just-prove-it.com